


## NE-382-1.7 -SRA , -SRA-NC Micro Pipet System 1.7 Assembly and Use


Remove needle from the syringe.




Install Repeating Adapter onto the syringe. Follow the instructions that came with the product.

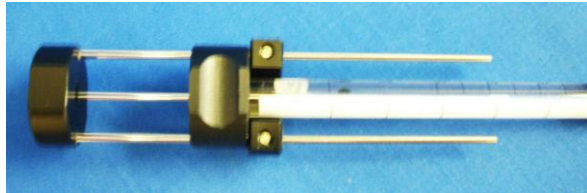

Insert the Bushing into the base of the Syringe in place of the needle. It will slip easily into position.

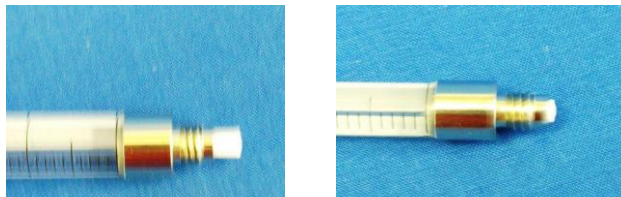


Thread the 2-piece Capillary Adapter onto the Syringe. See "How to assemble and use the Capillary Adapter NE-382-CAOR" that is provided on the next page.

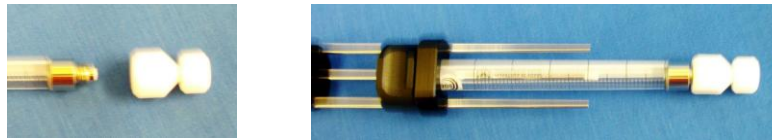

To insert the Glass "Needle" see "How to assemble and use the Capillary Adapter NE-382-CAOR" provided on the next page.




The Glass "Needles" are fragile-use caution and always use protective eyewear.

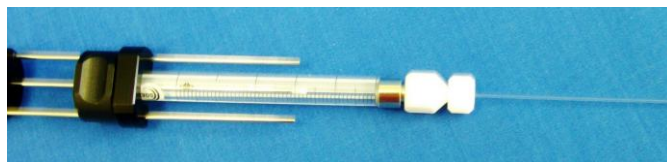


A: insert the 1.7mm capillary into the Capillary Holder.  
B: Place "needle" Guide over capillary

To transfer the sample, gently insert the "Needle" to the bottom of the capillary and slowly load the sample as you slowly lift the needle tip. Any air bubbles can be removed from the sample by spinning in a centrifuge.

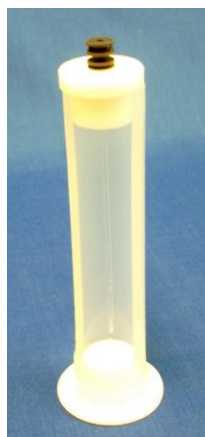

A

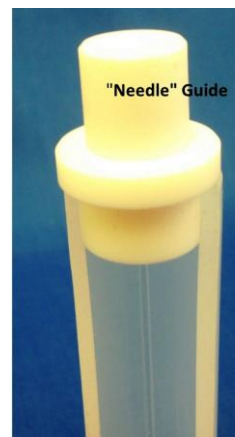

B